

**DO NOT** use this equipment unless an instructor has instructed you in its safe use/operation and has given permission.



Safety glasses may be required at times in work areas.



Long and loose hair must be contained or constrained.



Appropriate footwear with closed-toe must be worn.



Long-sleeve or protective clothing is required.



Rings and jewellery must not be worn.



Protective gloves may be required.

## PRE-OPERATIONAL SAFETY CHECKS

1. Check workspaces and walkways to ensure no slip/trip-hazards are present.
2. Check that all equipment components are in position and are operational.
3. Ensure you are familiar with the operation of the equipment.
4. Keep table and work area clear of all tools and debris.
5. Faulty equipment must not be used. Immediately report suspect equipment.
6. Be aware of any other students or personnel in the immediate vicinity and ensure the area is clear before using this equipment.
7. Familiarize yourself with all mechanical operations, switches and controls, including potential chemicals that may be used.
8. If you have any questions or concerns, ask the laboratory technician or instructor first.

## OPERATIONAL PROCEDURES

### **NO FOOD OR DRINK PERMITTED NEAR EXPERIMENTS.**

1. Turn on power switch in rear of equipment.
2. Fill up both containers with water.
3. Turn on water heater.
4. Adjust dial as needed.
5. Ensure water has reached boiling.
6. Take measurement readings.
7. Dispose of water afterwards.

## HOUSEKEEPING

1. Turn off all equipment and dispose of water on completion. Leave equipment and working area in a safe, clean and tidy state.

## POTENTIAL HAZARDS

- Steam burn on skin
- Electrical shock
- Manual handling
- Boiling water

Date of last review \_\_\_\_\_

Signature \_\_\_\_\_